



Terahertz Crystals

GaSe • ZnTe

SEMICONDUCTOR TERAHERTZ CRYSTALS

ZnTe

ZnTe (Zinc Telluride) crystals with $\langle 110 \rangle$ orientation are used for THz generation by optical rectification process. Optical rectification is a difference frequency generation in media with large second order susceptibility. For femtosecond laser pulses which have large bandwidth the frequency components interact with each other and their difference produce bandwidth from 0 to several THz.

Detection of the THz pulse occurs via free-space electro-optic detection in another $\langle 110 \rangle$ oriented ZnTe crystal. The THz

pulse and the visible pulse are propagated collinearly through the ZnTe crystal. The THz pulse induces a birefringence in ZnTe crystal which is read out by a linearly polarized visible pulse. When both the visible pulse and the THz pulse are in the crystal at the same time, the visible polarization will be rotated by the THz pulse. Using a $\lambda/4$ waveplate and a beamsplitting polarizer together with a set of balanced photodiodes, it is possible to map THz pulse amplitude by monitoring the visible pulse polarization rotation after the ZnTe crystal

at a variety of delay times with respect to the THz pulse. The ability to read out the full electric field, both amplitude and delay, is one of the attractive features of time-domain THz spectroscopy.

ZnTe are also used for IR optical components substrates and vacuum deposition.

NOTE: ZnTe crystal contains micro bubbles and they are visible in projection of illuminated crystal. However this does not affect the THz generation. We do not accept complains on presence of bubbles in crystal.



ZnTe, $\langle 110 \rangle$ CUT

Catalogue number	Size, mm	Thickness, mm	Holder, mm	Price, EUR
ZnTe-100H	10×10	0.1	Ø25.4	890
ZnTe-200H	10×10	0.2	Ø25.4	730
ZnTe-500H	10×10	0.5	Ø25.4	620
ZnTe-1000H	10×10	1.0	Ø25.4	510
ZnTe-2000H	10×10	2.0	Ø25.4	750

GaSe

GaSe (Gallium Selenide) crystals used for THz generation shows a large bandwidth of up to 41 THz. GaSe is a negative uniaxial layered semiconductor with a hexagonal structure of 62 m point group and a direct bandgap of 2.2 eV at 300 K. GaSe crystal features high damage threshold, large nonlinear optical coefficient (54 pm/V), suitable transparent

range, and low absorption coefficient, which make it an alternative solution for broadband mid infrared electromagnetic waves generation. Due to broadband THz generation and detection using a sub-20 fs laser source, GaSe emitter-detector system performance is considered to achieve comparable or even better results than using thin ZnTe crystals. In order to achieve

frequency selective THz wave generation and detection system, GaSe crystals of appropriate thickness should be used.

NOTE: because of material structure it is possible to cleave GaSe crystal along (001) plane only. Another disadvantage is softness and fragility of GaSe.



GaSe crystal mounted in Ø25.4 mm holder

GaSe, Z-CUT

Catalogue number	Clear aperture, mm	Thickness, µm	Holder, mm	Price, EUR
GaSe-10H1	Ø7	10	Ø25.4	1950
GaSe-30H1	Ø7	30	Ø25.4	1625
GaSe-100H1	Ø7	100	Ø25.4	1475
GaSe-500H1	Ø7	500	Ø25.4	1460
GaSe-1000H1	Ø7	1000	Ø25.4	1635
GaSe-2000H1	Ø7	2000	Ø25.4	1810

Please note that from now all standard GaSe crystals are provided mounted into Ø25.4 mm ring holders. Crystals could be mounted into Ø40 mm holders under your request.